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Measurement Facilities in the Infrared-Optics Laboratory

Infrared Radiation

Radiative transfer is important in many technical applications such as building physics, thermal insulation in the aerospace industry and combustion processes. ZAE Bayern has different measuring devices for determining the energy emitted by a surface and the heat transfer through a material.

Devices

Reflectance and transmittance are measured with an integrating sphere (Fig.1).

The emittance at high temperatures is determined by comparing the energy emitted by the sample and by a black body at the same temperature.

Two Fourier-Transformation-Infrared (FTIR) spectrometers and a diffraction spectrometer cover the spectral range between 0.25 μm and 40 μm , which includes the ultraviolet, visible and infrared regions.

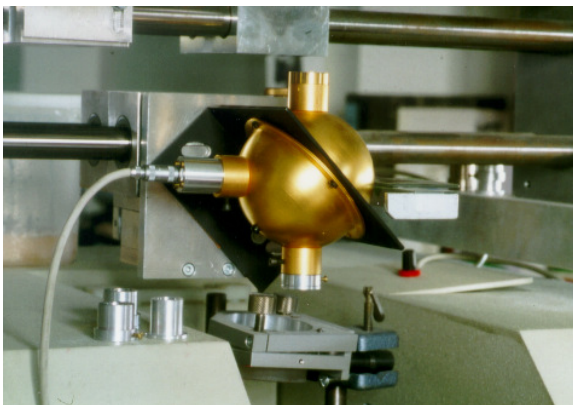
Fig. 1: FTIR spectrometer with integrating sphere.

Measurement Facilities

The devices at ZAE Bayern can measure:

- directional-hemispherical transmittance T_{dh} ,
- directional-directional transmittance T_{dd} ,
- directional-hemispherical reflectance R_{dh} ,
- directional-directional reflectance R_{dd} ,
- emittance ϵ .

The following wavelength ranges and temperatures are covered:

spectral values at room temperature		
value	wavelength	sample temperature
T_{dh}, R_{dh}, ϵ T_{dd}, R_{dd}	0.25 ... 500 μm	20 $^{\circ}\text{C}$

spectral values in dependence on temperature and emission angle		
value	wavelength	sample temperature
ϵ	2 ... 18 μm	200 ... 800 $^{\circ}\text{C}$

The actual spectral region in which measurements are carried out depends on the intended application and operation temperature.

Evaluation

Microscopic values, such as absorption and scattering, can be determined from the above-mentioned macroscopic values T_{dh} and R_{dh} with the aid of a special solution derived from the equation of radiative transfer. Absorption and scattering are characteristic for each material examined.

Furthermore, being able to calculate the mass-specific effective extinction e^* from the above enables the radiative thermal conductivity to be determined, which, in turn, enables the presentation of the radiative thermal transfer through materials. Mass-specific absorption a and scattering s^* can also be determined.

Contact:

Dr. J. Manara
Tel.: ++49-931/70564-46
e-mail: manara@zae.uni-wuerzburg.de
Dott.-Ing. M. Arduini-Schuster
Tel.: ++49-931/70564-17
e-mail: carla@zae.uni-wuerzburg.de
Fax: ++49-931/70564-60
<http://www.zae-bayern.de>