Permeations-Messtand

Motivation

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Gas Permeation System

Background

The permeability of foils and coatings is an important parameter in the fields of packaging, vacuum insulation and corrosion prevention. Knowing a material's selective permeability for different gases is a crucial aspect in energy-saving gas separation. Permeability is also a decisive characteristic for optimization in the fields of building physics, filter technology and porous electrodes. The permeameter system at ZAE Bayern can determine permeability for each of the above-mentioned applications within a range of approx. 15 decimal powers.

Measuring principle

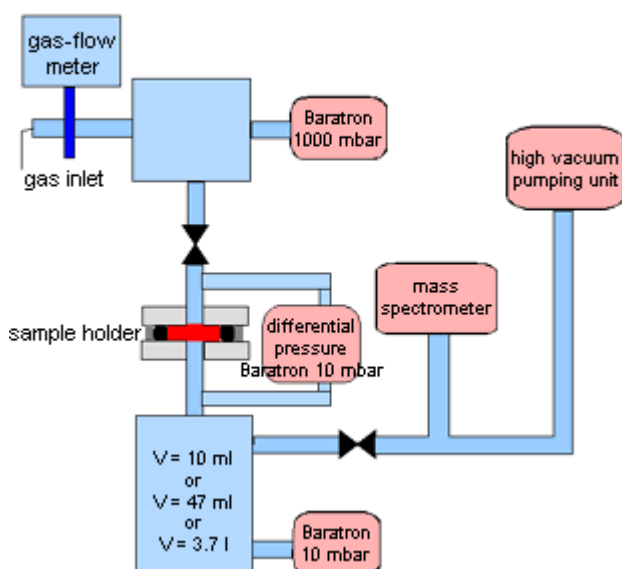
Fig. 1: Permeameter system – experimental set-up

There are two methods of carrying out measurements with the permeameter system (Fig. 1). On the one hand, the system enables the time-resolved determination of the integral via the gas-particle flow through a sample, whereby the pressure on the back of the sample is detected by a membrane pressure transducer after the measuring gas has been admitted to the front of the sample. Steady-state permeability measurements are also possible by determining the gas flow and pressure drop over the sample. This method is particularly useful for porous materials or materials with high permeabilities. On the other hand, the permeameter system can also differentially measure the gas-particle flow through the sample, whereby some of the atoms or molecules which penetrate the sample after the measuring gas has been admitted to the front of the sample are detected on the back of the sample by a mass spectrometer. The mass spectrometer provides a signal proportional

to the gas-particle flow. To calculate the gas-particle flow on an absolute scale, the system has to be calibrated with the aid of integral measurement. The mass spectrometer enables specific gas types to be detected, for example, when a mixture of gases is admitted to the sample. All the variables can be determined as a function of the average gas pressure between 10 mbar and 1000 mbar.

Specifications

Measurable variables: Permeability, solubility, adsorption behaviour

Measurement range:

$$5 - 10^{-8} \frac{\text{cm}^3(\text{STP}) \cdot \text{cm}}{\text{cm}^2 \cdot \text{s} \cdot \text{bar}} \cdot \frac{d(\text{cm})}{A(\text{cm}^2) \cdot \Delta p(\text{bar})}$$

whereby d = sample thickness,
 A = penetrated area,
 Δp = pressure gradient over the sample.

Temperature: 25 - 80 °C

Atmosphere: 5 - 1000 hPa

Measuring gases: He, Ar, N₂, N₂/O₂, CO₂, air

Sample sizes:

- Foils: $\varnothing = 40$ mm,
maximum thickness = 1 mm
- Discs: $\varnothing = 9 - 20$ mm,
maximum thickness = 3 mm
- Porous media: $\varnothing = 9 - 20$ mm,
maximum thickness = 3 mm

Contact:

Dr. G. Reichenauer
Tel.: ++49-931/70564-28
Fax: ++49-931/70564-60
E-mail: reichenauer@zae.uni-wuerzburg.de
<http://www.zae-bayern.de>

Address:

ZAE Bayern
Am Hubland
97074 Würzburg
Germany